ImageQuant TL 8.1 安裝及啟用流程

注意事項

一、此 IQTL 軟體為十人網路授權版為了不影響其他使用者權益,請分析完成 後必須離開程式,以利其他使用者登錄使用 二、若無法登錄即為使用人數已滿,請耐心稍後再行登錄

1.請打開 IQTL 資料匣



2.點選資料匣内 setup 檔案

| 2稱 | 修改日期 | 類型 | 大小 |
|---|--------------------|-------|--------|
| 📙 iqtl installation | 2017/8/31 下午 02:47 | 檔案資料夾 | |
| iqtlsecurity installation | 2017/8/31 下午 02:47 | 檔案資料夾 | |
| license server | 2017/8/31 下午 02:47 | 檔案資料夾 | |
| AutoRun | 2017/8/17 下午 04:01 | 安裝資訊 | 1 KB |
| Install | 2017/8/17 下午 04:01 | 園示 | 6 KB |
| IQTL 8.1 8.1.15 web installation instruct | 2017/8/17 下午 04:01 | 文字文件 | 2 KB |
| 🗿 Setup | 2017/8/17 下午 04:01 | 應用程式 | 138 KB |

3.點選 Install ImageQuant TL 方框, 即可安裝 ImageQuant TL 軟體 ℤ IQTL Software Installation

| mageQuant" TL | Install ImageQuant TL Getting Statted Readme User manual | Installed |
|---|---|---------------|
| noge Analyses Software vr.0 () () () () () () () () () () () () () (| Install IQTLSecuriTy | Not Installed |
| | Readme User manual | 1 |
| | Install License Server | Not Installed |
| imagination at work | License Activation Homepage | |
| ageQuant is a trademark of GE althcare Companies. Adobe and obe Reader are trademarks of Adobe stems Inc. | Computer Ethernet Address: Unknown | |
| y use of this software is subject to GE althcare Standard Software End-User ense Agreement for Life Sciences ftware Products | Install Adobe® Reader® | Not Installed |
| Healthcare Life Sciences AB, a neral Electric Company. | | |
| 3E Healthcare Life Sciences AB, a Seneral Electric Company. 9 2008 General Electric Company – All | | |

4.點擊 ImageQuant TL icon



5.點選 1D gel analysis

| IQ Control (| Centre Centre |
|--------------|--|
| Imag | geQuant TL |
| | <u>1D gel analysis</u> Analyze a 1D electrophoresis gel image |
| 9 | Analysis Toolbox Analyze an image using area and profile-based tools |
| | Colony Counting Count colonies or detect 2D electrophoresis spots |
| | <u>Array analysis</u> Analyze a dot/slot blot, microtiter plate or macroarray |
| ? | Online Help Look up queries in the comprehensive reference guide |
| (egg | Also contained in this installation: |
| | / <u>FluorSep</u> I <u>Q Tools</u> |

6.選擇要分析的圖檔

| 搜尋位置(I): | 2channel.d | ir | • | ÷E | * | |
|---|---|---------------------------------------|-----------------|---------------|----------|------------|
| 点面 媒體櫃 電腦 | 1d_Tiers Array_Blot Array_micro Colony_Pet Colony_Spo | otitre ri ots | | | | |
| 網路 | 檔案名稱(N): 檔案類型(T): | 選擇要進行分析之 All ImageQuantTL image fi | 圖檔 es (*.tif | ;*.tiff;*.gel | • | Open 取消 |
| Analyse this image E 1D for analysis of mar | e using: | ✓ Preview | | | | |
| electrophoresis ge | á. | Preview not avails | ıble. | | | |

7.選擇 "I am using a Floating E-License"

| I am using a Node Locked E-license I am using a Floating E-License | | | |
|--|---------------------------------|-------|--|
| I am using a Floating E-License | 🔍 I am using a Node Locked E-li | tense | |
| . | I am using a Floating E-Licens | e | |
| | | | |
| | | | |
| | | | |

8.在空格內輸入 IP address: 120.126.98.236 後點選 Finish 即可使用

| ImageQuantTL License Setup Wizard | $\sim c$ | X | | | |
|---|-----------------------------|----------------------|--|--|--|
| License Server: 120.126.49.16 前入IP: 120.12 | 26.98.236 | | | | |
| If you do not have a license file, please collect one using t available on the GE Healthcare E-Licensing Webpage www.elicensing.amershambiosciences.com/gt/we | the Software Licensing Syst | em <mark>tool</mark> | | | |
| Please see the IQTL Getting Started guide for instructions on how to set up a license server. MAC address for this computer: 002618d2beed | | | | | |
| < Back Finish | Help | Quit | | | |